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
High-performance, Low-Toxicity Quantum Dots

— Next-Generation Fluorescent Probes for Biomarkers —

Cadmium- and selenium-free quantum dots with outstanding stability compared to organic dyes.

● What are Quantum Dots?

- Quantum dots are semiconductor nanoparticles approximately 2–10 nm in size, and their emission wavelength varies depending on the particle diameter.
- They exhibit exceptionally high photostability and are resistant to photobleaching even under prolonged illumination.
- Although they have been commercialized for displays, conventional quantum dots contain cadmium or selenium, requiring handling as hazardous materials.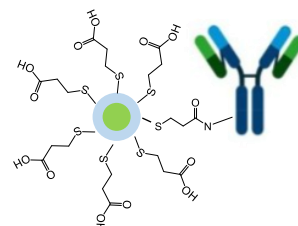

● Key Features of Our Development

◆ Non-toxic


These water-dispersible quantum dots are based on indium phosphide (InP) and achieve high performance without containing any toxic elements such as cadmium or selenium. They do not require disposal as hazardous materials, making handling significantly easier. In addition, cellular cytotoxicity testing confirmed $IC_{50} > 100 \mu\text{g/mL}$, indicating a high level of safety for biological applications.

◆ High Reproducibility


Quantum dots have traditionally been regarded as having low lot-to-lot reproducibility, but through precise control of the particle surface state, we have achieved highly reproducible reactivity with biological materials. The carboxyl groups introduced on the particle surface allow stable conjugation of antibodies and other biomolecules. Furthermore, pre-conjugated secondary-antibody quantum dots are available, enabling users to begin evaluation immediately.

◆ Easy Fluorescence Imaging

Quantum dots exhibit exceptional photostability, showing minimal photobleaching even under strong excitation—far superior to conventional organic dyes. Their absorption spectra extend continuously toward shorter wavelengths, allowing flexible selection of excitation wavelengths without strict consideration of the Stokes shift. *Note: Care should be taken regarding autofluorescence when imaging biological samples.*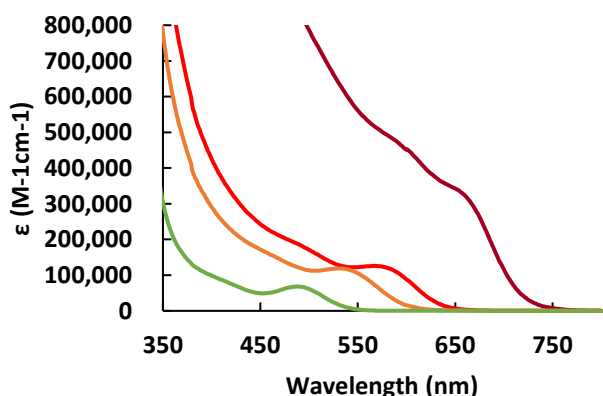

Molar absorption coefficient per particle

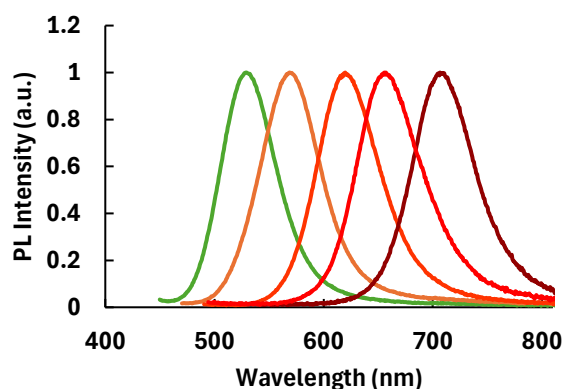Wide PL spectral range

● Example of Application

- Fluorescence imaging, in vitro diagnostics, flow cytometry, and other research applications


NIPPON CHEMICAL INDUSTRIAL

Organic Materials R&D Department
 Tel. +81-3-3636-8089
 Email: tomo.sakanoue@nippon-chem.co.jp




● Photobleaching Comparison in Confocal Laser Microscopy

We compared the photobleaching resistance of Alexa Fluor 488 with that of 710-nm-emitting quantum dots. Both probes were conjugated to anti-EGFR antibodies and used to stain A431 cells, followed by continuous imaging under 488-nm laser excitation. Alexa Fluor 488 showed rapid photobleaching even at low excitation power, whereas the quantum dots exhibited almost no photobleaching, even when the excitation intensity was increased. These results suggest that quantum dots enable more diverse and flexible imaging strategies—such as strong excitation, long-exposure imaging, and signal accumulation—fluorescent dyes. that are difficult to achieve with conventional dyes.

Alexa Fluor488-anti EGFR Antibody
Ex.=488 nm, Em.=525-545 nm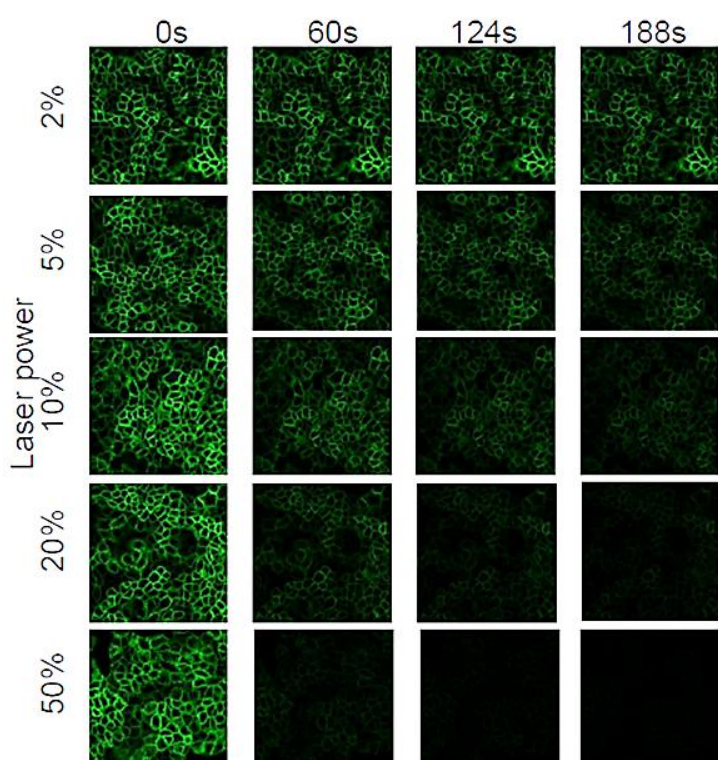

InP-QDs (710 nm) -anti EGFR Antibody
Ex.=488 nm, Em.=695-715 nm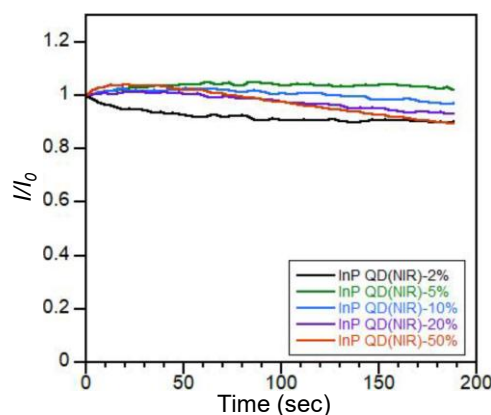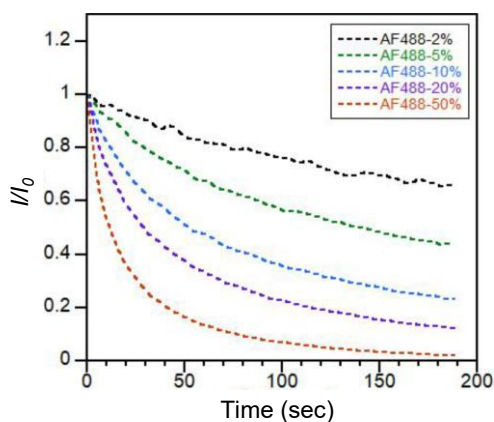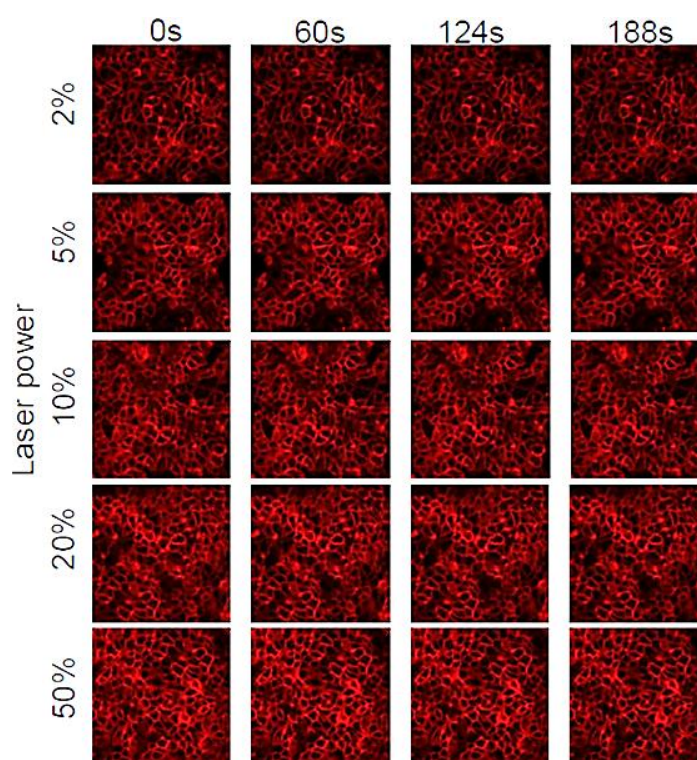

Courtesy of Ms. Yuna Sano, Univ. of Tokyo (Evident FV4000)

- These quantum dots are being developed in collaboration with Assoc. Prof. S. Ohta at Univ. of Tokyo.